

## Handling & Cleaning of Trays - Ambiset & BioFix

### Handling Instructions:

- Always hold the tray by the metal side handles while transporting.
- Do not carry the tray while the lid is open.
- Ensure the tray is placed on a flat, stable surface before opening.
- To open, unlock the latches and gently lift the lid until it reaches its stopping point.
- When closing, ensure that both the latches are fully shut and securely locked.

### Cleaning Instructions (as per use):

- Place the tray on a flat surface, such as a counter, or in a sink
- Open the lid and run warm water over the tray for several seconds.
- Gently pat the tray dry.
- Allow to air dry or initiate the “Tray Dry” cycle on the Stabilizer
- If using the “Tray Dry” cycle, close and lock the tray lid, place the tray inside the Stabilizer, and run the cycle. The Tray Dry cycle runs for 20 minutes.

### Storage Instructions:

- Store the trays in a clean environment.
- Ensure the tray is stored flat, with the mesh side facing down.
- Avoid stacking any items on top of the trays during storage.

### Cleaning the Insert Plates:

- Clean the insert plates with water or an ethanol spray as required.
- Dry the plates gently using lab wipes.

